



Job Aid

Document G_10_J_2_A Author: C Allen Authorised: W Porau 07/09/22 Review: 07/09/24

How to use a ZeoMed Biohazard Spill Kit

The purpose of a spill kit is to contain & clean up laboratory spills of blood, & body fluids.

1. Remove contents of ZeoMed Biohazard spill kit.



2. Put on plastic apron, facemask with eyeshield and gloves.




3. Open one of the clinical waste bags Press firmly on the base pushing mins outer edges of the bag to stand it upright.



4. Open ZeoMed absorbent powder sachet and apply over the spill- allow 1-2 to absorb excess moisture and odours



5. Using the scoop and scraper provided, pick up the substance; place into the yellow contaminated waste bag with the scoop and scraper. Seal the bag securely with the tie provided. Open the second clinical waste bag, place the first sealed clinical waste bag inside the second waste bag.

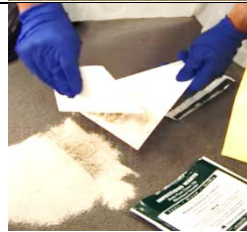

6. Clean the area with the detergent provided using one of the cleaning cloths. Put all cleaning cloths used into second clinical waste bag.



7. Remove apron, facemask & gloves. Place into the second clinical waste bag. Seal the bag securely with a tie provided. Dispose of waste in a contaminated waste bin.



8. Wash and dry hands thoroughly.

9. Notify the Lab Manager of the incident and complete an incident report form supplied by your employer. **Reorder Spill kit refill.**

Reference video https://www.youtube.com/watch?v=FD_sEU7vpMA

